Mikroskop · Microscope · Microscopio

Science ETD-201

Art. No. 5806200

DE Schnellstartanleitung

EN Quickstart guide

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Deutschland

<http://www.bresser.de>

For information regarding liability claims or service requests, please refer to chapters "Warranty" and "Service" in this documentation. We ask for your understanding, that any requests or submissions that have been addressed to the manufacturer's address can not be processed.

Errors reserved - technical specifications subject to change.

Copyright: 2016Bresser GmbH

All rights reserved.

It is not allowed to reproduce this documentation or parts of it in any kind (e.g. photocopying, printing, etc) without written a permission of the manufacturer. It is also prohibited to use or transmit this documentation with electronic systems (e.g. image file, website, etc).

The descriptions and brand names of the corresponding firms used in this documentation are generally protected under trademark, brand and/or patent law for Germany, for the European Union and/or other countries.

2 About this Instruction Manual

NOTICE

This Quick Start guide is not intended to replace the more detailed user manual.

Please read the safety instructions and the operating instructions carefully before use.

3 Intended Use

- This device is not intended for use by individuals (including children) with limited physical, sensory or mental capabilities or those lacking in experience and/or knowledge, unless they are supervised by an individual responsible for their safety or have received instructions from them regarding the use of the device.
- This product is intended only for private use.
- It was developed for the magnified display of things in nature.
- The device is intended only for indoor use.

4 General safety instructions

⚠ DANGER

Danger of an electric shock!

This device contains electronic components that operate via a power source (power supply and/or batteries). In case of any improper use of this device, there is a risk of an electric shock. An electric shock can cause severe injury or even death. Therefore please read the safety instructions below to avoid an explosion.

- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Always pull on the plug to separate the device from the power supply. Never pull on the cable.
- Before operating, check the device, cables and connections for damage.
- Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Only use the device in complete dry environment and do not touch it with wet or moist parts of your body.

DANGER

Danger of choking!

In case of any improper use of this device, there is a risk of choking, especially for children. Therefore please read the safety instructions below.

- Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard!

DANGER

Danger of explosion!

In case of any improper use of this device, there is a risk of an explosion. Therefore please read the safety instructions below to avoid an explosion.

- Do not expose the device to high temperatures. Use only the supplied power adapter. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

CAUTION

Danger of injury!

This device contains components and/or accessories that can cause minor to severe injuries in case of any improper use. Therefore please read the safety instructions below to avoid any bodily injury.

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

CAUTION

Fire hazard!

In case of any improper use of this device, there is a risk of fire. Therefore please read the safety instructions below to avoid the initiation of burning.

- Never cover the ventilation slots or cooling fins of the device while using it or as long as it has not sufficiently cooled down!

NOTICE**Risk of property damage!**

In case of any improper use of this device and/or its accessories, there is a risk of property damage. Therefore only use the device according to the safety instructions below.

- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose this device to higher temperatures and protect it from water and high humidity.
- Protect the device from severe shocks!
- For this device only use accessories and spare parts that comply with the technical information.

5 Product images

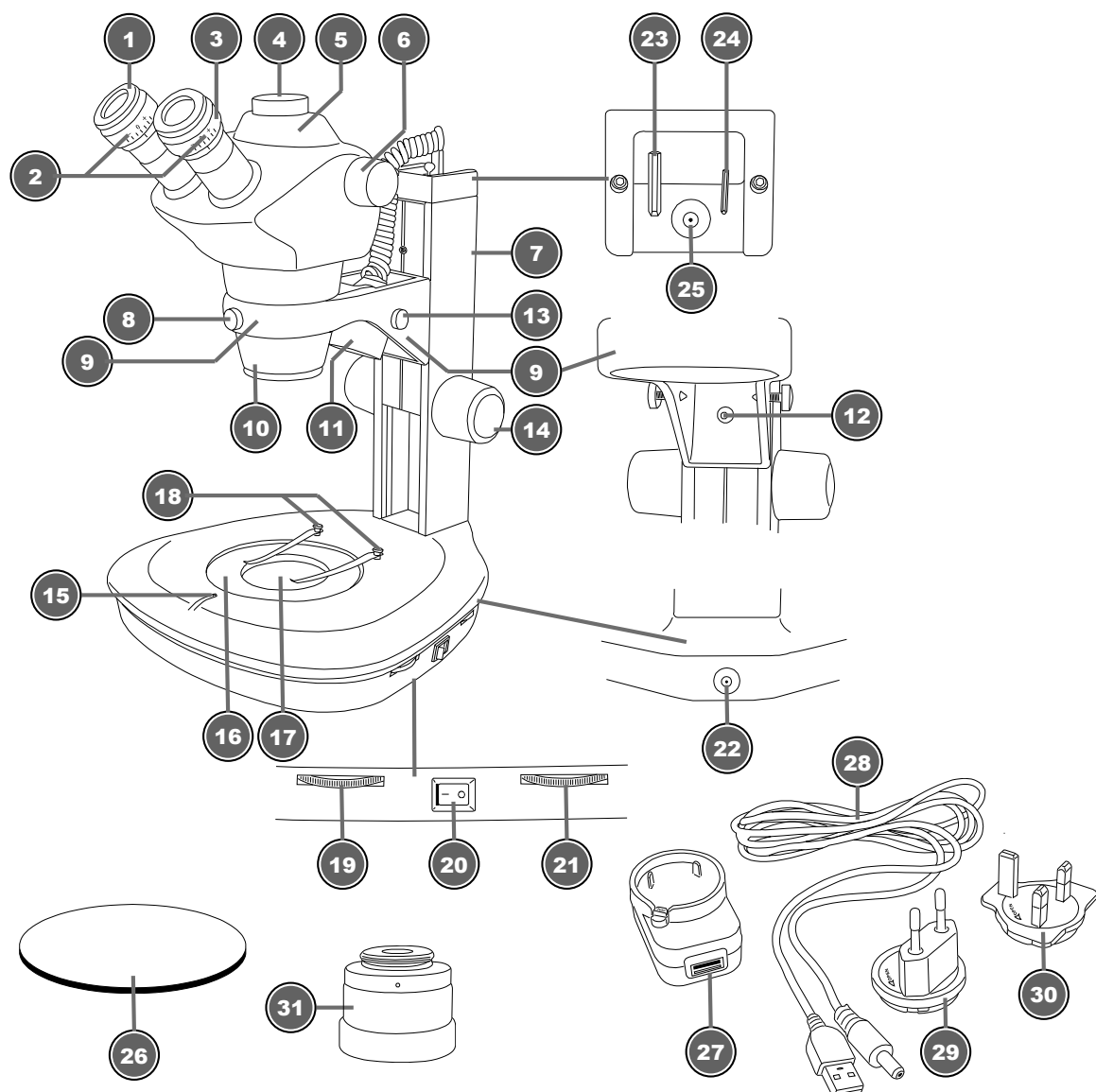
Illustration 1: Product images

6 Parts overview

1	Eyecup
2	Eyepiece
3	Diopter adjustment ring
4	Trinocular head
5	Microscope head
6	Zoom knob
7	Microscope stand
8	Fixing screw objective tube
9	Microscope head bracket
10	Objective tube with safety glass
11	Incident light illumination
12	Fixing screw microscope head bracket
13	Fixing screw incident light illumination
14	Focusing knob
15	Locking screw object plate
16	Object plate (glass)
17	Transmitted light illumination
18	Object fixing clamps
19	Transmitted light dimmer
20	On/Off switch
21	Incident light dimmer
22	DC connection socket for power cable
23	Allen wrench, large
24	Allen wrench, small
25	DC connection socket for incident light illumination with rubber cover
26	Two-colour object plate
27	Universal USB power adapter
28	USB power cable with coaxial/barrel connector
29	Power adapter attachment (EU)
30	Power adapter attachment (UK)
31	Photo adapter (C-mount)

7 Site location

Choose a suitable location before assembly and commissioning.

Place the device on a stable and flat surface, free from vibrations.

8 Installing the USB power adapter

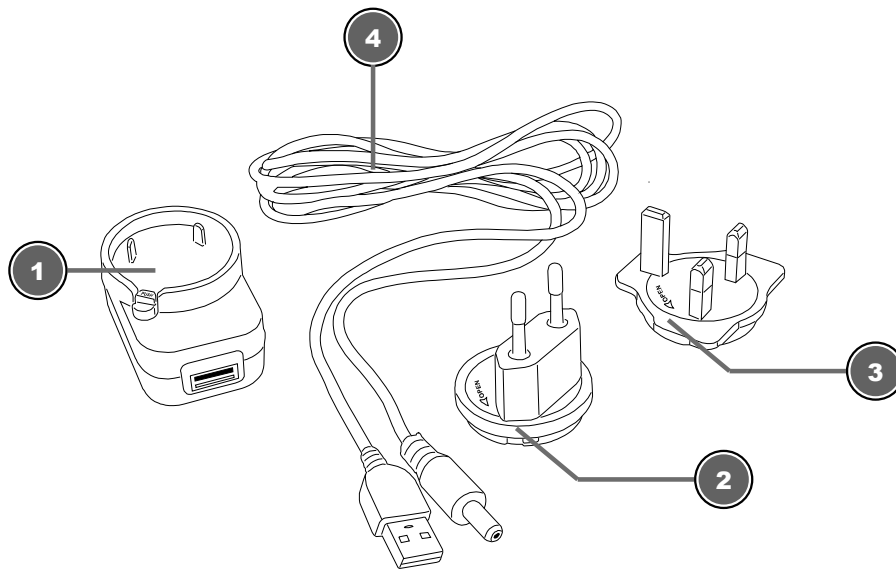
Illustration 2: USB power adapter with exchangeable attachment

1 USB power adapter	2 Adapter attachment (EU)
3 Adapter attachment (UK)	4 USB power cable with coaxial/barrel connector

1. Place suitable adapter attachment with its notches on the pins of the USB power adapter.
2. Turn adapter attachment in LOCK direction until it is locked.
3. Put the USB plug of the power cable into the USB port of the power adapter.
4. To remove the adapter attachment, push the PUSH button and turn the attachment in OPEN direction.

9 Turning on/off the device

1. Move On/Off switch to position 'I' to turn on the device.
2. Move On/Off switch to 'O' position to turn off the device.

10 Adjusting the lighting

This device is equipped with two lighting units with LED light sources which can be dimmed independently to each other.

To control the lighting units proceed as follows:

1. Turn on the device.
2. Turn/Move the dimmer control of the desired light source to adjust its brightness.

To adjust the angle of the light source, proceed as follows:

1. Loosen on of the locking screw for the incident light illumination.
2. Move the lamp holder to the desired position and re-tighten the locking screw by hand.

11 Cleaning and maintenance

- Protect the device against dust and moisture.
- Avoid putting fingerprints and equal contaminations on any optical surfaces.
- Make sure the residual moisture is fully eliminated before use.

- Disconnect the device from the power supply by pulling the power plug when it is not used or in case of longer interruption of operation and before starting any work on maintenance and cleaning.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.
- Do not use organic solvents (e.g. alcohol, ether, acetone, xylol or other dilutions) to clean lacquered parts or plastic parts!
- Disassemble optical elements (e.g. objective, eyepiece, etc) before cleaning.
- Blow away loose dust from the lens surfaces first.
- Use high quality lens cleaning towels or soft cloth and moisten it with a little bit of pure alcohol (available in a drug store). Wipe the lens surface with it.
- Pull a dust protection cover* over the microscope and store it in a closed container at a dry and mould-free place. Store the microscope and the accessories in the relevant containers when they are not used for a longer time. It is recommended to store also objectives and eyepieces in closed containers with desiccant.

(*depending on the model a dust protection cover and box can be part of the content)

NOTICE

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

12 Warranty

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

13 EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards.

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

e-mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd
Unit 1 starborough Farm,
Starborough Road, Nr Marsh Green,
Edenbridge, Kent TN8 5RB
Great Britain

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Folux B.V.