
DIGITALES MIKROSKOP DIGITAL MICROSCOPE 20X/80X/350X

- DE** Bedienungsanleitung
- EN** Operating instructions
- FR** Mode d'emploi
- NL** Handleiding
- IT** Istruzioni per l'uso
- ES** Instrucciones de uso
- RU** Руководство по эксплуатации

Lieferumfang · Contents · Contenu de la livraison · Inhoud · Accessori in dotazione
Artículos incluidos en la entrega · Material fornecido · Комплект поставки

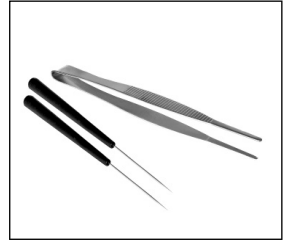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Name: Telefon:

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General Warning

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.
- This device contains electronic components which operate via a power source (power supply and/or batteries). Only use the device as described in the manual, or you run the risk of an electric shock.
- Do not expose the device to high temperatures. Use only the power adapter supplied. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or even an explosion.
- Never bend, pinch, pull or run over the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as they pose a choking hazard.
- Before cleaning the device, disconnect the device from the power supply by removing the plug.
- Position your device so that it can be disconnected from the power supply at any time. The wall socket you use should be located near the device and easily accessible, since the plug on the power cable serves as a disconnecting device for the power supply.
- Always pull on the plug to separate the device from the power supply. Never pull on the cable.
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to temperatures above 45°C.

Declaration of Conformity

Bresser GmbH has issued a 'Declaration of Conformity' in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request.

The Parts of your Microscope

1. Microscope body
2. Camera shutter release
3. Magnification adjustment knob
4. Microscope camera
5. Surface illumination
6. Camera mount
7. Microscope stage
8. Substage illumination
9. Slide clips
10. Slide
11. Focus adjustment knobs
12. Microscope stand foot
13. Metal spring for electrical contact
14. Guide bar for electrical contact
15. Buttons for adjusting the brightness (+/-)
16. On/off button for type of illumination
17. USB port on the microscope
18. Slide box with 5 permanent slide preparations, 5 empty slides and 10 cover slips
19. Set of instruments for microscope
20. Dust cover
21. Software CD
22. USB connection cable

1. Assembling the Digital Microscope

Before you can use your microscope for the first time, you must assemble it properly. First, remove all parts from the packaging. Place the microscope body (1) with the stand foot (12) on a table. Screw the microscope camera (4) into the camera mount (6) from above.

Important:

When inserting the camera, be sure that the metal spring (13) on the camera (4) leads into the guide bar (14) on the camera mount (6). Plug the USB connection cable (22) into the USB port on the microscope (17), inserting the end with the smaller connector.

2. Using the Software CD

You'll find a software CD with the driver and image-processing software packaged along with your new microscope. You can install the software on your computer. When you connect your microscope to your computer with the USB cable, you can look at the pictures on the computer monitor and save them using the 'CamApp' software. You can edit the pictures with the 'Photomizer' software. You'll learn how to properly install the software in the next sections.

Note:

In order to work with the Digital Microscope for the long term, we recommend that you always connect it to the same USB port.

2.1 Installation of the Software

Important: Before inserting the CD into your computer, first plug the USB cable into the USB port on your PC. Windows will recognize the new device and will indicate this in a notification window. Click on 'Cancel', since the driver for the device will be installed automatically later.

1. Save all current data and close any open programs.
2. Now insert the CD-ROM that came with your microscope into the CD/DVD drive of your computer. The installation menu starts automatically. If it does not, go to the Windows explorer and select the CD/DVD drive (most of the time, it's the 'D' drive, but it can have another letter). From there, start the file 'menu.exe' by double-clicking with the left mouse button.

2.2 Driver installation

To install the driver software, click on 'Install Driver' on the menu with the mouse cursor. Then follow the installation programme instructions. During the software installation the correct driver for your operating system will be installed automatically. No manual input is needed. Once the computer restarts, the driver is installed. In rare cases the device may not be recognised by your computer. If this occurs, you need to install the driver manually from the CD. Select the menu item 'Repair Program' during the installation of the driver. After the driver has been installed, restart the computer. If this fails, please refer to the troubleshooting chapter that follows.

2.3 Installation of the Image Editing Software Photomizer

The image editing software 'Photomizer' is located on the software CD. You can edit your pictures here.

Note: The 'Photomizer' software requires 'Microsoft .NET Framework 3.5'. This must be installed first, if it has not been installed on your operating system already.

You can directly download this here:

<http://www.microsoft.com/downloads/en/details.aspx?FamilyID=333325fd-ae52-4e35-b531-508d977d32a6>

Please note that the manufacturer cannot guarantee the accessibility of the site or that it is free from viruses. Likewise, they cannot guarantee the functionality of the software.

1. To install the software, click once on the menu item 'Install Software' with the left mouse button.
2. A selection window appears in which you can select the preferred language. Select the preferred language and confirm with 'OK'.
3. In the welcome window, please click on 'Continue'.
4. In the next window, you will be asked for the 'Destination Folder'. Here, just click on 'Next'.
5. Click on 'Continue' in the 'User-defined Installation' window.
6. The options 'Create Desktop Shortcut' or 'Link in the Quick Launch Bar' set up program shortcuts on your desktop and taskbar.
7. Now, the window with the setup status will appear. A progress bar will inform you about the ongoing installation. This process can take a few minutes.
8. The window 'Photomizer is being installed' appears. Click on 'Finish'. The installation ends.

2.4 Troubleshooting: The device isn't recognised by the computer (USB).

1. This occurs in rare cases with some OEM Windows Vista or XP versions. The device is not defective. Updating the notebook or PC USB driver usually fixes this problem.
2. If, after installation of the driver, the computer tries to operate the device on another USB port, the driver may need to be re-installed from the CD. Select the menu item 'Repair Program' during the installation of the driver. After the driver has been installed, restart the computer.

3. If not, proceed as below. (The maker cannot accept any liability whatsoever in these cases¹). Delete ghost devices. These are devices not actually currently connected to your computer. The reason ghost devices appear is that Windows generates a system settings entry for each USB device (e.g. USB drive) on each port. Windows then seeks that device each time it's started. This slows down USB device recognition and Windows booting and causes nonrecognition of USB ports. Cleaning up the system environment and deleting devices that are no longer in use helps here. To do so, right click 'Workplace' ('Computer' in Vista) and open the 'Properties' menu. Then click the 'Advanced' tab ('Advanced System Settings' in Vista) and then 'Environment Variables'. Choose the 'New' option. Enter the following under 'Variable Name': **devmgr_show_nonpresent_devices**. Under 'Variable Value' enter '1'. Confirm your input with 'OK' and then reboot the computer. Change to the device manager. In the 'View' menu activate the 'Show Devices Option'. The ghost devices previously not shown will now be displayed in light grey. Check the various categories including USB and memory size. Only delete entries for devices you no longer use.¹

¹Source: Microsoft Knowledge Base

<http://support.microsoft.com/kb/315539/en>

3. Working with the Digital Microscope

3.1 Setting the illumination

You do not need any extra electrical cord for the light, because the Digital Microscope gets the electricity for the light via the USB cable. The Digital Microscope offers you three illumination options:

- a) surface illumination (light from above),
- b) substage illumination (light from below) or
- c) surface and substage illumination
(light from above and below at the same time).

Their use depends on the object you would like to observe (see section 3.4). If you press the on/off button (16) once, the surface illumination (5) will be turned on. If you press the button a second time, the illumination switches to substage (8). If you then press the button a third time, both surface and substage illumination are engaged. Another press of the button turns the lighting completely off again. Choosing the right lighting depends on what you would like to look at (see section 3.4).

3.2 Live Observation

If you press the camera shutter release (2) on your microscope, the 'CamApp' window opens. If the program 'CamApp' does not start automatically, switch to 'All Programs'. You can start the 'CamApp' program manually under 'USB2.0 PC Camera'.

A live image (in general, blurry) is displayed on the monitor. Focus the live image by turning the setting wheels for the image sharpness (11).

3.3 Choosing the Magnification Settings

You should begin your first observation with the smallest magnification (20X). This makes adjusting the image much easier, because the higher the magnification, the more carefully the object of observation must be positioned (with the transmitted-light microscope, for instance, a slide would be the object of observation). Therefore, turn the magnification adjustment knob (3) to "20X". Then, adjust the microscope stage (7) to its lowest position using the focus adjustment knobs (11).

3.4 Observing an Object

When observing objects through a microscope, you must distinguish between transparent and opaque objects. Transparent objects are generally completely flat or liquid (e.g. insect wings, a drop from a puddle, etc.); opaque objects are generally larger and not so flat (e.g. smaller insects, plant pollen, rocks, coins, etc.).

3.4.1 Observing Transparent Objects

Your Digital Microscope comes with permanent slide preparations. These are slides on which transparent objects have already been mounted and prepared for observation with a microscope. Set one of these permanent slide preparations onto the microscope stage (7). Push the slide underneath both slide clips on the microscope stage, so that nothing can move about during the observation. Then, turn the light to substage illumination (8) by pressing the on/off button for the light (16) as described in section 3.1. Next, adjust the permanent slide so that the observation object is clearly visible on the monitor, in the preview area of the camera software. Now, slowly turn the focus adjustment knob (11) left or right until the image is in focus. By pushing multiple times on the buttons for light brightness (+/-) (15), you can set the illumination so that the fine details of the observation object appear clearly on the monitor. If you would like to switch to a higher magnification, turn the magnification adjustment knob (3) to the left, to the position 80X or 350X.

Important:

Each time you adjust the magnification, you must also use the magnification adjustment knob (3) to change the focus.

3.4.2 Observing Opaque Objects

If you would like to observe opaque objects with this microscope, the object must be lit from above. Therefore, use the on/off button for the illumination (16) to set the lighting to surface illumination (5), as described in section 3.1. Place the observation object on the small, round disk in the middle of the microscope stage. Using the focus adjustment knob (11), focus the image. It will also appear in focus on the monitor. You might have to move the observation object around a little bit so that it is displayed properly on the monitor. By pushing multiple times on the buttons for light brightness (+/-) (15), you can set the illumination so that the fine details of the observation object appear clearly on the monitor. The Digital Microscope allows you to observe objects with a height of up to 6 mm.

Tip:

If you are observing objects that are only slightly transparent (e.g. minerals, grains of sand, etc.), you can improve the picture quality by using surface and substage illumination at the same time (section 3.1.). The brightness for both lighting types is set with the appropriate keys (+/-) (15).

Important:

If you turn on the illumination on the Digital Microscope in a very bright room or near a window on a sunny day, the image on the monitor can appear much too bright. This is because the camera is very sensitive. Too much light overpowers the fine details of the observation object. In this case, you must completely shut off the light on the Digital Microscope, so that all the details are visible on the monitor.

3.5 Taking Pictures

Using the camera shutter (2), you can take a picture and save it as a JPG file.

1. Press the camera shutter (2) or select 'Still image capture' in the menu 'Capture(R)' in the 'CamApp' program.
2. The 'Photo' window appears on the screen with a picture.
3. To save the image, click 'File' and 'Save' or 'Save As'.

3.6 Making Movies

The 'CamApp' program allows you to make movies with the microscope and save them as AVI files.

1. By choosing 'File' from the menu, you can specify a folder to save your AVI files to under 'Set Capture File Folder'. For example, 'Own Files' -> 'Own Videos'.
2. You can prepare to film with 'Capture(R)' and start filming with 'Start Capture' from the menu.
3. Under the 'Capture' menu, you can end your filming with 'Stop Capture'.
4. You can watch your film on your PC monitor using a playback program for multimedia files.

4. Starting and Using the 'Photomizer' Software

1. Start the 'Photomizer' software on your computer.
2. In the program window, click 'Open Camera'.
3. If you have more than one device connected to your PC, you can choose the desired device in the next step. If you only have one connected device, you can skip this step.
4. Now, you should be able to see the image from the camera on your monitor. On the microscope, bring the image into focus by turning the focus adjustment knobs (11).
5. Click 'Capture' in order to capture an image you would like to save. This will be shown on the right in the menu bar.
6. Select this image with a single mouse click, then click on 'Transfer Image'.
7. You will now leave the image-capture stage, and go back to the program window.
8. Now click on 'File' at the top of the screen and then on 'Save File' in order to save the image.

Tip:

If you have questions or problems with the 'Photomizer' software, you can always go to the program's menu bar, and click on '?' and then on 'Open Help'. For questions or problems, you can look on the website www.photomizer.net. Please ask your parents before using the Internet.

5. Accessories

Slides: Place the specimen you wish to observe on a slide. After you have covered it with some water or gum media, you can place a cover slip on top.

Cover slips: Use the cover slips to cover specimens that you have placed on a slide in order to protect them while you observe them with the microscope.

Tweezers: These are a type of pincers that you can use to pick up small objects.

Dissecting needle: A dissecting needle can be used for many things. You can use it to scratch the surface of a specimen, to fix objects in place or to stir liquids.

6. Care and maintenance

The microscope is a high-quality optical appliance. Therefore you should make sure that dust and moisture do not come into contact with your microscope. Avoid putting fingerprints on any optical surfaces.

If dust or dirt has gotten onto your microscope or the accessories, you should first remove the dirt with a soft brush. Then clean the soiled area with a soft, lint-free cloth. To remove fingerprints from the optical surfaces, it is best to use a lint-free, soft cloth, to which you have applied some alcohol.

7. Technical Data

LEDs: 5 pieces

System requirements: PC with Intel Pentium IV processor or higher; Operating System: Windows XP, Windows Vista or Windows 7; at least 1024 MB RAM memory; at least 500 MB free hard disk space; free USB 2.0 port; DirectX; Service Pack 3 for Windows XP, Service Pack 2 for Windows Vista; .NET Framework 3.5; Internet Access

Warranty and warranty term extension

The warranty term is two years from the date of purchase. Please keep your proof of purchase. Register at www.bresser.de/warranty and fill out a brief questionnaire to get your warranty term extended to **five years**. Registration must be completed within three months of purchase (date of receipt) to validate the warranty. If you register thereafter, the warranty term will not be extended. If you have problems with your device, please contact our customer service first. Do not send any products without consulting us first by telephone. Many problems with your device can be solved over the phone. If the problem occurred after the warranty ended or it is not covered by our warranty terms, you will receive a free estimate of repair costs.

Service Hotline: +49 (0) 2872 - 80 74-210

Important for any returns:

Please make sure to return the device carefully packed in the original packaging in order to prevent damage during transport. Also, please enclose your receipt for the device (or a copy) and a description of the defect. This warranty does not imply any restriction of your statutory rights.

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are trademarks of the National Geographic Society,
used under license. All rights reserved.