# BIOLUX NV

Art. No. 96-19150

- Ⓧ DE Bedienungsanleitung
- Ⓧ GB Operating Instructions
- Ⓧ FR Mode d'emploi
- Ⓧ NL Handleiding
- Ⓧ IT Istruzioni per l'uso
- Ⓧ ES Instrucciones de uso
- Ⓧ PT Instruções de utilização
- Ⓧ FI Käyttöohje
- Ⓧ SE Bruksanvisning
- Ⓧ SI Navodila za uporabo
- Ⓧ HU Használati utasítás
- Ⓧ SK Návod na použitie
- Ⓧ CZ Návod k použití
- Ⓧ PL Instrukcja obsługi
- Ⓧ RO Instrucțiuni de folosire

Fig. 1





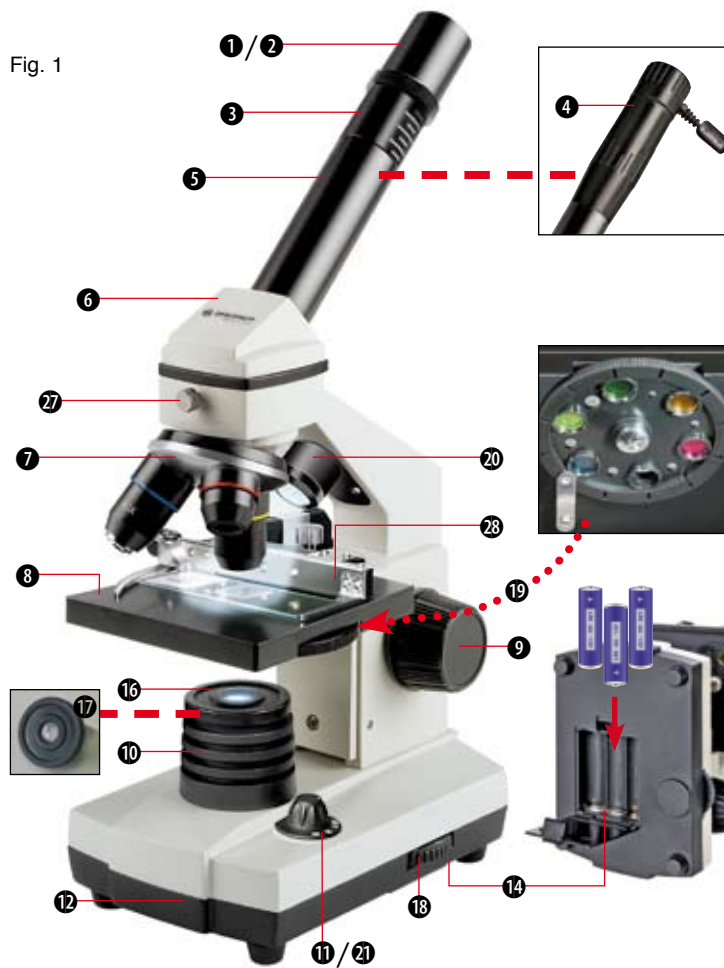
Fig. 2
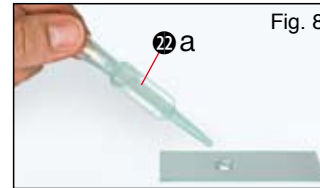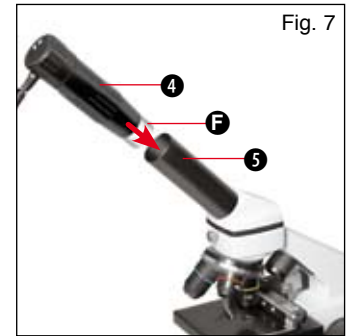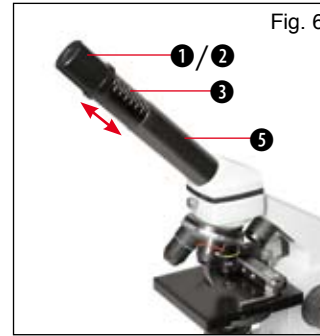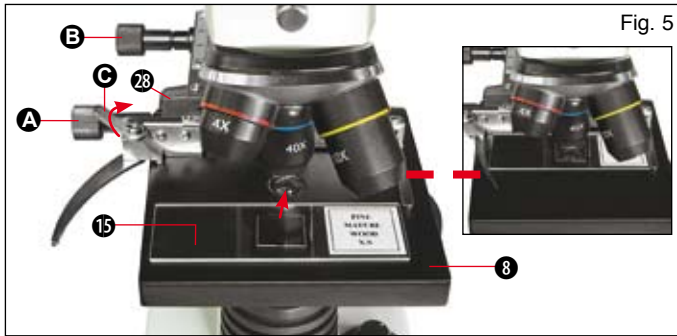
<b>DE</b>	<b>Bedienungsanleitung .....</b>	<b>5</b>
<b>GB</b>	<b>Operating Instructions.....</b>	<b>17</b>
<b>FR</b>	<b>Mode d'emploi.....</b>	<b>28</b>
<b>NL</b>	<b>Handleiding.....</b>	<b>40</b>
<b>IT</b>	<b>Istruzioni per l'uso.....</b>	<b>51</b>
<b>ES</b>	<b>Instrucciones de uso .....</b>	<b>63</b>
<b>PT</b>	<b>Instruções de utilização .....</b>	<b>75</b>
<b>FI</b>	<b>Käyttöohje .....</b>	<b>87</b>
<b>SE</b>	<b>Bruksanvisning .....</b>	<b>97</b>
<b>SI</b>	<b>Navodila za uporabo.....</b>	<b>108</b>
<b>HU</b>	<b>Használati utasítás.....</b>	<b>119</b>
<b>SK</b>	<b>Návod na použitie .....</b>	<b>130</b>
<b>CZ</b>	<b>Návod k použití.....</b>	<b>141</b>
<b>PL</b>	<b>Instrukcja obsługi.....</b>	<b>152</b>
<b>RO</b>	<b>Instrucțiuni de folosire.....</b>	<b>164</b>

---

## Allgemeine Informationen

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e-Mail:   service@bresser.de

Rhede, 2013-11-01

A handwritten signature in blue ink, appearing to read 'Helmut Ebbert'.

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## General Information

### About this Instruction Manual

These operating instructions are to be considered a component of the device. Please read the safety instructions and the operating instructions carefully before use. Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

#### **DANGER!**

You will find this symbol in front of every section of text which deals with the risk of severe injury or even death in the event of improper use.

#### **CAREFUL!**

You will find this symbol before every section of text that deals with the risk of minor to severe injuries resulting from improper use.

#### **NOTE!**

You will find this symbol in front of every section of text which deals with the risk of damage to property or the environment.

### Intended Use

This product is intended only for private use. It was developed for the magnified display of things in nature.

## General Warning

### The device is intended only for indoor use.

#### **DANGER!**

Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

This product contains small parts which could be swallowed by children. There is a RISK OF CHOKING.

This device contains electronic components which operate via a power source (batteries). Children should only use the device under adult supervision. Only use the device as described in the manual; otherwise, you run the risk of an electric shock.

Do not expose the device to high temperatures. Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.

Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.

Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat.

Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

Children should only use the device under adult supervision. Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard.

#### **CAREFUL!**

Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after use. In the event of contact with the eyes or mouth, rinse thoroughly with water. In the event of pain, contact a doctor immediately and take the substances with you.

Before cleaning the device, disconnect it from the power supply by removing the plug or batteries.

## **!** NOTE!

Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. The batteries should be removed from the unit if it has not been used for a long time.

Never recharge normal, non-rechargeable batteries. This could lead to explosion during the charging process.

Do not disassemble the device. In the event of a defect, please contact the Service Centre in your country (see the warranty card).

## Operating instructions

### Parts overview (Fig. 1+2):

- ❶ 5x WF Eyepiece
- ❷ 16x WF Eyepiece
- ❸ Barlow lens
- ❹ MicrOcular (PC-Ocular)
- ❺ Eyepiece holder
- ❻ Microscope head
- ❼ Objective-revolver
- ❽ Microscope stage
- ❾ Focus wheel
- ❿ LED lighting (transmitted light)
- ⓫ Illumination ON/OFF
- ⓬ Microscope base
- ⓭ Photomizer software
- ⓮ Battery compartment
- ⓯ 5 Slides, 10 Covering glasses and 5 preparations in a plastic box
- ⓰ matted lens
- ⓱ Condenser lens
- ⓲ Dimmer
- ⓳ Colour filter disc
- ⓴ LED lighting (direct light)
- ⓵ Direct light / transmitted light switch
- ⓶ Microscope tools:
  - a) Pipette; b) Tweezers
- ⓷ Prawn breeding plant
- ⓸ MicroCut
- ⓹ Specimens:
  - a) yeast; b) Gum media (specimen

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- ⓺ Carrying case
- ⓻ Locking screw
- ⓼ Mechanical desk

### 1. General/Location

First you must make sure that your microscope is on a stable and solid surface.

### 2. Electric LED lighting with dimmer

Check before use whether the light switch (Fig. 1, 21) is set to "off".

The microscope has two lighting units. Lighting can be of three types. Set the switch (Fig 1, 21) to „II“ to light the specimen from above (reflected light) or „I“ to light it from below (direct light). Use setting „III“ to have the specimen simultaneously subjected to direct and transmitted light. The transmitted light unit (Fig 1, 10) is used for transparent specimens (those on glass slides). To view solid non-transparent specimens use the direct light unit (Fig 1, 20). Use of both forms of lighting simultaneously is only recommended for semi-transparent specimens. This operating mode is not recommended for direct light specimens on slides as it may cause reflection on the slide.

The batteries supplied should be inserted into the battery compartment located in the base plate before use. (Fig. 1. 14)

Remove the battery compartment cover by pressing lightly on the cover clip. Insert the batteries into the holder.

Important! Pay attention to the polarity as marked on the battery holder terminals.

As your device has continuously controllable lighting (dimmer) optimal illumination of the object to be viewed is guaranteed.

### 3. Colour filter disc

The colour filter under the microscope table (Fig. 1, 19) aids in viewing very bright and transparent objects. Just select the right colour for the specimen in question. The components of colourless/transparent objects (e.g. starch particles, single-cell -specimens) can thus be better recognised.

### 4. Interchangeable illumination lenses

The illumination of your microscope contains two lenses (Fig. 2, 16+17). These are – according to the object being viewed – to be attached to the LED illumination for best results. The ground glass lens Fig. 1, 16) is already mounted on the lighting unit. To change the lenses, simply screw them on and off. Just turn the upper lighting unit (Fig. 1, 10).

An overview of purposes:

**Matted lens (Fig. 2, 16)** to be used for

- views with the PC eyepiece (Fig. 2, 4)

- viewing of extremely small items with eyepiece (Fig. 2, 1+2) and Barlow lens (Fig. 2, 3)

**Condenser lens (Fig. 6, 17)** to be used for

- viewing of standard items with eyepieces (Fig. 2, 1+2) and Barlow lens (Fig. 2, 3)

### 5. Attach mechanical desk

A mechanical desk is included along with your microscope (located in the case, under the CD). To install this on the microscope, proceed as follows:

1. Place the mechanical desk (Fig. 4, 28) on the microscope stage (Fig. 4, 8) in such a way that the attachment screws (Fig. 4, D) are positioned over the middle threaded hole (Fig. 4, E).
2. Tighten the attachment screws securely.

### 6. Microscope setup

The microscope's eyepiece (Fig 1, 6) will now be prepared for the first observation.

First, loosen the screw (Fig. 1, X27) and rotate the eyepiece into a convenient position.

Begin every observation with the lowest magnification.

Place the microscope's table (Fig. 1, 8) with the focus knob (Fig 1, 9) into the lowest position and rotate the objective revolver (Fig. 1, 7) until it locks on the lowest magnification (4x).

#### **NOTE:**

First, drive the Microscope's table (Fig. 1, 8) to its lowest position before changing the objective in order to prevent damage.

Insert the 5x eyepiece (Fig. 3, 1) in the Barlow lens (Fig. 3, 3).

Take care, that the Barlow lens is inserted completely in the monocular head (Fig. 3, 5).

### 7. Observation

After you have set up the microscope with the corresponding illumination, the following principles are important:

Begin each observation with a simple observation at lowest magnification, so that the centre and position of the object to be viewed is in focus.

The higher the magnification the more light is required for good picture quality.

Place a permanent slide culture (Fig. 5, 15) directly under the microscope lens on the plate (Fig. 5, 8) and clamp it on the cross-table (Fig. 5, 28). To do so push the lever (Fig. 5, C) aside. The specimen to be examined must be directly over the lighting. If not turn the two knurled screws (Fig. 5, A+B) on the mechanical plate.

**TIP:**

On the mechanical plate (Fig 5, 28) there are two knurled screws (Figs 5, A+B). They are used to precisely position the specimen laterally (Fig 5, A) and vertically (Fig 5, B).

Look through the eyepiece (Fig. 1, 1/2) and turn carefully the focusing wheel (Fig. 1, 9) until you can see a sharp picture.

Now you can get a higher magnification, while you pull out slowly the Barlow lens (Fig. 6, 3) of the monoculare barrel (Fig. 6, 5). With the nearly entirely pulled out Barlow lens the magnification is raised to 2x.

For still higher magnification you can put the 16x eyepiece (Fig. 2, 2) into the objective revolver (Fig. 1, 7) and set on higher position (10x / 40x).

**TIP:**

Depending on the preparation higher magnifications do not always lead to better pictures.

With changing magnification (eyepiece or objective lens changes, pulling out of the Barlow lens) the sharpness of the image must be newly defined by turning the focusing wheel (Fig. 1, 9).

**NOTE:**

Please be very careful when doing this. When you move the mechanical plate upwards too fast the objective lens and the slide can touch and become damaged.

## 8. Viewed Object – condition and preparation

### 8.1. Condition

With the Barlow lens nearly fully extended magnification can be doubled. Both transparent and non-transparent specimens can be examined with this microscope, which is a direct as well as transmitted light model. If opaque specimens are examined - such as small animals, plant parts, tissue, stone and so on - the light is reflected from the specimen through the lens and eyepiece, where it is magnified, to the eye (reflected light principle, switch position I). If opaque specimens are examined the light from below goes through the specimen, lens and eyepiece to the eye and is magnified en route (direct light principle, switch position II). Many small water organisms, plant parts and the finest animal components have natural transparent characteristics others need to be prepared accordingly. Preparation

can take many forms such as the soaking in appropriate substances (medium) to make the object opaque. Objects can also be finely sliced (hand cut, micro-cut). The following sections explain these procedures in detail.

### 8.2. Creation of thin preparation cuts

Specimens should be sliced as thin as possible, as stated before. A little wax or paraffin is needed to achieve the best results. A candle can be used for the purpose. The wax is put in a bowl and heated over a flame. The specimen is then dipped several times in the liquid wax. The wax is finally allowed to harden. Use a MicroCut (Fig 2, 24) or knife/scalpel (carefully) to make very thin slices of the object in its wax casing.

**DANGER!**

Be extremely careful when using the knives/scalpels or the MicroCut. There is an increased risk of injury due to the sharp edges!

These slices are then laid on a glass slide and covered with another.

### 8.3. Creation of an own preparation

Put the object to be observed on a glass slide and add a drop of distilled water using a pipette (Fig. 8, 22a) onto the object (Fig. 8).

Set a cover glass (available in most well stocked hobby shops) perpendicularly at the edge of the water drop, so that the water runs along the cover glass edge (Fig. 8). Now lower the cover glass slowly over the water drop.

**TIP:**

The gum medium supplied (Fig 2, 25b) is used to make permanent slide cultures. Add it instead of distilled water. The gum medium hardens so that the specimen is permanently affixed to its slide.

## 9. Experiments

If you have made yourself familiar with the microscope already, you can accomplish the following experiments and observe the results under your microscope.

### 9.1. Newspaper print

**Objects:**

1. A small piece of paper from a newspaper with parts of a picture and some letters
2. A similar piece of paper from an illustrated magazine

Use your microscope at the lowest magnification and use the preparation of the daily paper. The letters seen are broken up,

because the newspaper is printed on raw, inferior paper. Letters of the magazines appear smoother and more complete. The picture of the daily paper consists of many small points, which appear somewhat dirty. The pixels (raster points) of the magazine appear sharply.

### 9.2. Textile fibres

**Objects and accessories:**

1. Threads of different textiles: Cotton, line, wool, silk, Celanese, nylon etc..
2. Two needles

Each thread is put on a glass slide and frayed with the help of the two needles. The threads are dampened and covered with a cover glass. The microscope is adjusted to a low magnification. Cotton staples are of vegetable origin and look under the microscope like a flat, turned volume. The fibres are thicker and rounder at the edges than in the centre. Cotton staples consist primary of long, collapsed tubes. Linen fibres are also vegetable origin; they are round and run in straight lines direction. The fibres shine like silk and exhibit countless swellings at the fibre pipe. Silk is animal origin and consists of solid fibres of smaller diameter contrary to the hollow vegetable fibres. Each fibre is smooth and even moderate and has the appearance of a small glass rod. Wool fibres are also animal origin; the surface consists of

overlapping cases, which appear broken and wavy. If it is possible, compare wool fibres of different weaving mills. Consider thereby the different appearance of the fibres. Experts can determine from it the country of origin of wool. Celanese is like the name says, artificially manufactured by a long chemical process. All fibres show hard, dark lines on the smooth, shining surface. The fibres ripple themselves/crinkle after drying in the same condition. Observe the thing in common and differences.

### 9.3. Salt water prawns

**Accessories:**

1. Prawn eggs (Fig 2, 25d)
2. Sea salt (Fig 2, 25c)
3. Prawn breeding plant (Fig 2, 23)
4. Yeast (Fig 2, 25a)

**CAREFUL!**

Eggs and prawns are not fit for human consumption.

### 9.3.1 The lifecycle of the saltwater prawn

The saltwater prawn or *artimia salina* to scientists has an unusual and interesting lifecycle. The female's eggs are hatched without any male shrimp having to fertilise them. The resultant baby prawns are all female. Under unusual circumstances such

as when a swamp is drained the eggs may produce male prawns. These males fertilise the female's eggs, resulting in a specific type of eggs. These are called winter eggs and have a thick shell as protection. They're pretty rugged and can survive the swamp or lake drying out causing the death of the entire prawn population for up to a decade in a form of hibernation. The eggs hatch once again as soon as the correct ambient conditions have been obtained. The eggs supplied (Fig 2, 25d) are of this type.

### 9.3.2. Hatching of the salt water prawn

To hatch the prawn it is essential to first have a saline solution suited to the prawn's needs. Fill half a litre of rain- or fresh-water in a container. Let it stand for about thirty hours. As water evaporates over time it's a good idea to have a second container of such water left standing for thirty-six hours. Once stood for this length of time pour half of the sea salt supplied into one of the containers (Fig 2, 25c) and stir until it has dissolved. Then pour some of it into the prawn breeding plant (Fig 2, 23). Add a few eggs and close the lid. Put it somewhere with plenty of light but not in the direct sun. The temperature should be approximately 25° C. The prawns will hatch in two or three days at this temperature. Should any water evaporate during this time replace it from the second container.

### 9.3.3. The saltwater prawn under the microscope

What comes out of the egg is known as a nauplius larva. Use the pipette (Fig 2, 22a) to put some of them on a slide for examination. They will move in the solution using their hair like limbs. Remove a few daily from the container for examination under the microscope. If you do so and save the pictures made with the MicrOcular you will then have a seamless record of the prawn's lifecycle. You can remove the upper lid of the prawn bleeding plant and put the whole thing under the microscope. The larvae will mature in six to ten weeks depending on ambient temperature. You will soon have bred an entire generation of saltwater prawns that constantly reproduce.

### 9.3.4. Feeding your saltwater prawns

To keep them alive saltwater prawns must be fed occasionally. This must be done carefully as overfeeding causes the water to stagnate and poison the prawns. Feeding is best done with dry powdered yeast (Fig 2, 25a). Give them a little every other day. If the water darkens this signifies it is stagnating. If so remove the prawns and put them in a fresh saline solution.

## 10. MicrOcular setup

### NOTE:

The MicrOcular only works without the Barlow lens supplied. Magnification setting is changed by using the MicrOcular and must be reset by re-focussing.

Remove the Barlow lens (Fig 6, 3) and eyepiece currently in use from the eyepiece holder (Fig 6, 5) and replace them with the MicrOcular (PC-Ocular) (Fig 7, 4) and reducer lens (Fig 7, F) as in illustration 7 in the holders (Fig 7, 5)

### NOTE:

Please do not yet connect MicrOcular and PC. Please go through the items below in sequence.

## 11. Installation and use of the software

### 11.1. Software and installation information

A software CD is included with your microscope. The software and drivers on this CD must be installed on your computer in order to use your MicrOcular. Once you connect the MicOcular to your computer after

installation, you can view pictures on and save them to your computer using Photomizer. To install the software and driver correctly, simply follow the installation steps.

## 11.2. Software installation

1. **Important:** Before inserting the CD, first plug the USB cable into the USB port on your PC. Windows will now recognize the new device, and will indicate this in a notification window. Now please click on "Cancel".
2. Now insert the CD-ROM that came in your package into the CD/DVD drive of your computer. The installation menu starts automatically. If it does not, go to the Windows explorer and select the CD/DVD drive (most of the time, it's the "D" drive, but it can have another letter). From there, start the file "autorun.exe" by double-clicking with the left mouse button.

### 11.2.1. Driver installation

To install the driver software, click on the menu point "Install Driver" with the mouse cursor. Then follow the installation program instructions. During the software installation the correct driver for your operating system will be installed automatically. No manual input is needed. In rare cases the device may not be recognised by your computer. As a rule you need then only install the driver manually from the CD. If this

fails please refer to the troubleshooting chapter that follows.

### 11.2.2. Installation of the Image Editing Software Photomizer Pro (Bresser Edition)

The image editing software "Photomizer" is located on the software CD. You can edit your pictures here.

1. To install the software, click once on the menu point "Install PHOTOMIZER" with the left mouse button.
2. The Photomizer Software requires Microsoft .NET Framework 4.0, which it will install if it is not already on your system. If it is already on your system, you may skip to step 5.
3. In the welcome window, please accept the Microsoft license agreement, and then click "Install". The installation can take a few minutes.
4. As soon as everything is installed, click "Finish".
5. Now you will be presented with a choice, in which you can choose your language. Make your selection and confirm it by clicking „OK“.
6. When you see "Welcome", click on „Next“.
7. In the next window, you will be asked for the „Destination Folder“. Here, just click on „Next“.

8. Now, the window with the Setup-Status will appear – here, a progress bar will inform you about the ongoing installation. This process can take a few minutes.
9. The window "Photomizer is being installed" appears. Click on "Finish". The installation ends.

#### **TIP:**

To use the MikrOkular long-term we recommend it always be connected to the same USB port.

## 12. Using the MicrOcular

### 12.1. Preparation

1. Slide a specimen under your microscope and focus on it.
2. Remove the eyepiece and Barlow lens from the eyepiece support and the dust cap from your MikrOkular and install same in the eyepiece supports instead of the Barlow lens.
3. Start your PC if you haven't yet and connect your MikrOkular to the USB port of your computer.

### 12.2. Showing and saving MikrOkular images on your PC

1. Start the Photomizer Software.
2. Click „Open camera“

3. In the event that you have connected more than one unit, you can choose the desired unit in the subsequent selection. Click on "USB 2.0 Webcam". If only one device is connected, this step is omitted.
4. The camera image should now be visible on your screen. Focus the microscope image.
5. Click „Capture“ to record an image. It will then be shown on the right.
6. Click on it to select it and then click „Transfer image“.
7. This takes you to the Photomizer software.
8. File - Save as

### 12.3. The Photomizer Pro Software (Bresser Edition)

If you need help in the „Photomizer“ programme click „?“ and then „Open help“. If you need further assistance please visit the maker's home page at [www.photomizer.net/bresser](http://www.photomizer.net/bresser)

## Care and maintenance

Separate the device from the power supply before cleaning (remove batteries)!

Only use a dry cloth to clean the exterior of the device.

#### **NOTE:**

Do not use any cleaning fluid to avoid damaging the electronics.

Clean the lenses (eyepiece and/or lens) with a soft and lint-free cloth only (e.g. microfibre).

#### **NOTE:**

Do not apply excess pressure to the cloth so as to avoid scratching the lenses.

To remove more stubborn dirt, moisten the cleaning cloth with an eyeglass-cleaning solution and wipe the lenses gently.

Protect the device from dust and moisture. Store it in the supplied bag or transportation packaging. The batteries should be removed from the unit if it has not been used for a long time.

## Troubleshooting

### **Error**

No picture visible

### **Solution**

switch light on  
put condenser lens  
in place  
adjust focus

Picture flickers  
while viewing  
with PC eyepiece

if necessary,  
adjust resolution of the  
video graphics board

Software installation  
reports  
“not XP approved”

confirm by  
clicking “OK”

## Specifications

### **System requirements for PC eyepiece**

Minimum system requirements: PC with an Intel Pentium IV processor or higher; Windows XP with Service Pack 3\*, Windows Vista (32/64Bit) with Service Pack 2\* or Windows 7 (32/64Bit) with Service Pack 1\*; .NET Framework 4.0\*; min. 1024 MB RAM (64Bit = 2048 MB); min. 500 MB free hard drive space; free USB port; CD/DVD/BD drive.

\*available for free via Windows Update (Internet connection needed)

## Magnification table

<i>Eyep.</i>	<i>Objct.</i>	<i>Magn.</i>	<i>with Barlow I.</i>
5x	4x	20x	40x
5x	10x	50x	100x
5x	40x	200x	400x
16x	4x	64x	128x
16x	10x	160x	320x
16x	40x	640x	1280x

## Disposal

Dispose of the packaging materials properly, according to their type (paper, cardboard, etc). Contact your local waste disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household waste!

As per the Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

In accordance with the regulations concerning batteries and rechargeable batteries, disposing of them in the normal household waste is explicitly forbidden. Please make sure to dispose of your used batteries

as required by law – at a local collection point or in the retail market. Disposal in domestic waste violates the Battery Directive.

Batteries that contain toxins are marked with a sign and a chemical symbol. „Cd“ = cadmium, „Hg“ = mercury, „Pb“ = lead.

<sup>1</sup> battery contains cadmium

<sup>2</sup> battery contains mercury

<sup>3</sup> battery contains lead

## SERVICE AND GUARANTEE

This device is guaranteed for 5 years from the date of purchase. Please keep the receipt as proof of purchase. Our products are manufactured using the latest manufacturing techniques and undergo rigorous quality control.

### Service Centre

Please contact the service centre for any questions regarding the product or claims. Our trained staff will be happy to help. The service hotline is listed below. If it is necessary to make a guarantee claim (repair or exchange), our service staff will supply you with the return address.

Service Hotline\*: 00800 6343 7000

### Correction of defects and returns

We will repair any material or manufacturing defects free of charge during the guarantee period. Excluded from the guarantee are faults due to damage caused by the user, or improper use due to non-authorized third-party attempted repairs or similar. If it is necessary to repair or exchange a product, our service centre employees will supply you with the return address.

In the case of returns, please note the following:

- Make sure that the article is carefully packaged. If possible use the original packaging.
- Fill in the returns form and include it with a copy of your receipt and the product you are returning.

## Returns form

Transaction no. \*\*: .....

Product name: .....

Short description of defect: .....

.....

.....

Surname, first name: .....

Postcode / Town: .....

Street / House no.: .....

Telephone: .....

Date of purchase: .....

Signature: .....

\*free call / \*\*will be supplied to you by the service centre

## Declaration of Conformity

Product Type: Microscope

Product Name: BRESSER Biolux NV

Article No.: 96-19150

Meade Instruments Europe GmbH & Co. KG declares that the above-named product corresponds to the following directives and relevant standards in its conception and design as well in the version marketed by us.

**Directive:**  
EMC 2004/108/EC

ErP 2009/125/EC

**Harmonized Standards:**  
EN 55011:2007+A2:2007  
EN 61326-1:2006  
Regulation (EC) 278/2009

This declaration is no longer valid in the event of a change to the product which has not been coordinated with us.

Meade Instruments Europe GmbH & Co. KG  
Gutenbergstr. 2  
D-46414 Rhede  
Germany  
e-Mail: [service@bresser.de](mailto:service@bresser.de)

Rhede, 2013-11-01

A handwritten signature in blue ink, appearing to read 'Helmut Ebbert'.

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