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# Table of contents

<b>1</b>	<b>Imprint</b> .....	<b>18</b>
<b>2</b>	<b>Validity note</b> .....	<b>18</b>
<b>3</b>	<b>About this Instruction Manual</b> .....	<b>19</b>
<b>4</b>	<b>Intended use</b> .....	<b>19</b>
<b>5</b>	<b>General safety instructions</b> .....	<b>19</b>
<b>6</b>	<b>About this microscope</b> .....	<b>21</b>
<b>7</b>	<b>Features</b> .....	<b>21</b>
<b>8</b>	<b>Software Download</b> .....	<b>21</b>
<b>9</b>	<b>Choice of location</b> .....	<b>22</b>
<b>10</b>	<b>Parts overview and scope of delivery</b> .....	<b>22</b>
<b>11</b>	<b>Assembling the microscope</b> .....	<b>23</b>
<b>12</b>	<b>Power supply</b> .....	<b>23</b>
<b>13</b>	<b>Adjusting the lighting</b> .....	<b>23</b>
<b>14</b>	<b>Connecting to a PC</b> .....	<b>24</b>
<b>15</b>	<b>Adjust stand</b> .....	<b>24</b>
<b>16</b>	<b>Adjusting the zoom</b> .....	<b>25</b>
<b>17</b>	<b>Focusing the object</b> .....	<b>25</b>
<b>18</b>	<b>Recording</b> .....	<b>25</b>
<b>19</b>	<b>Technical data</b> .....	<b>26</b>

<b>20</b>	<b>CE declaration of conformity .....</b>	<b>26</b>
<b>21</b>	<b>UKCA declaration of conformity .....</b>	<b>26</b>
<b>22</b>	<b>Disposal.....</b>	<b>27</b>

---

# 1 Imprint

Bresser GmbH

Gutenbergstr. 2  
46414 Rhede  
Germany

[www.bresser.de](http://www.bresser.de)

For any warranty claims or service inquiries, please refer to the information on "Warranty" and "Service" in this documentation. We ask for your understanding that unsolicited returns cannot be processed.

Errors and technical changes excepted.

© 2023 Bresser GmbH

All rights reserved.

The reproduction of this documentation - even in extracts - in any form (e.g. photocopy, print, etc.) as well as the use and distribution by means of electronic systems (e.g. image file, website, etc.) without the prior written permission of the manufacturer is prohibited.

The designations and brand names of the respective companies used in this documentation are generally protected by trade, trademark and/or patent law in Germany, the European Union and/or other countries.

## 2 Validity note

This documentation is valid for the products with the following article numbers:

9652200 9652201 9652202

**Manual version:** 0324

**Manual designation:**

Manual\_9652200-L467811\_Digital-Microscope\_en-de-fr-it-es-nl-cs-sk\_BRESSER\_v032024a

Always provide information when requesting service.

---

### 3 About this Instruction Manual

---

#### NOTICE

**These operating instructions are to be considered a component of the device.**

Read the safety instructions and the instruction manual carefully before using this device.

Keep these instruction manual in a safe place for future reference. If the device is sold or passed on, the instruction manual must be passed on to any subsequent owner/user of the product.

---

### 4 Intended use

- This device is not intended for use by persons (including children) with impaired physical, sensory or mental abilities, or for lack of experience and/or knowledge, unless they are supervised by or have received instructions from a person responsible for their safety on how to use this device.
- This device is for personal use only.
- It was developed for an magnified representation of nature observations.
- The device is intended only for indoor use.

### 5 General safety instructions

---

#### DANGER

##### **Danger of suffocation!**

Improper use of this product may result in suffocation, especially for children. It is therefore imperative that you observe the following safety information.

---

- Keep packaging materials (plastic bags, rubber bands, etc.) away from children! There is a danger of choking!
-

- 
- This product contains small parts that can be swallowed by children! Choking hazard!
  - If small parts are swallowed, consult a doctor immediately!

## CAUTION

### Risk of body injury!

This device contains components and/or accessories that can cause minor or serious body injury if used improperly. It is therefore essential that you observe the following safety information in order to prevent body injuries.

- Sharp-edged and pointed tools are often used for working with this device. Therefore, keep this device and all accessories and tools in a place inaccessible to children. Danger of body injury!
- Chemicals and liquids do not belong in children's hands! Don't drink! Clean hands thoroughly under running water after use. In case of accidental contact with eyes or mouth, rinse with water. In case of complaints, consult a doctor immediately and show the substances.

## NOTICE

### Danger of material damage!

Improper handling may result in damage to the unit and/or accessories. Therefore, use the device only in accordance with the following safety information.

- Do not disassemble the device! In the event of a defect, please contact your dealer. They will contact the Service Center and can arrange the return of this device for repair if necessary.
- Do not expose the device to high temperatures and protect it from water and high humidity.
- Do not subject the device to excessive vibrations.

- 
- Only use accessories and spare parts for this device that comply with the technical specifications.

## 6 About this microscope

This microscope is only a digital microscope and therefore does not offer the possibility of visual observation.

It is equipped with an objective lens and a camera sensor for image recording. The digital images can be transferred directly to a Windows PC or Mac via the supplied USB cable, being viewed there and evaluated, optimized and finally saved using the supplied software\*.

\*Software available as download for Windows and Mac (see Software Download)

## 7 Features

- USB Digital microscope
- Bright ring light illumination from eight adjustable LEDs
- High quality zoom optics: 50x to 1000x magnification
- Working distance: 15mm (50x) to 40mm (1000x)
- Image format: BMP/JPG
- Image resolution: 640x480, 1280x720, 1920x1080\*
- Video format: AVI (30fps)
- PC interface: USB (5V)
- Incl. Windows PC software (download), stand

\*Interpolated

## 8 Software Download

### **HiView Software Download**

Required operating system: Windows (Windows 7 or later) or macOS (10.13 or later)

<http://www.bresser.de/download/9652200>

Alternate download:

<http://soft.hvscam.com>

## 9 Choice of location

Choose a suitable location before installation and commissioning.

Place the device on a stable, level and vibration-free surface.

## 10 Parts overview and scope of delivery

1 Digital microscope holder

2 Bracket for flexible alignment

---

3	Locking screw (holder)	4	Stand
5	Digital microscope	6	Dimmer button
7	Record button	8	Zoom wheel (50x to 1000x)
9	LED ring lighting	10	Objective
11	Protection cap	12	USB cable
13	USB-A connector		

### **Scope of delivery:**

Digital microscope (A) with USB cable, microscope stand (B)

## 11 Assembling the microscope

For the assembly proceed as follows:

Snap the microscope into the center of the clamp on the stand.

## 12 Power supply

Plug the USB-A connector of the USB cable into a free USB port of the PC. The power is supplied through the computer after it has been switched on. The LED ring lighting will light up.

## 13 Adjusting the lighting

The digital microscope is equipped with a LED ring lighting for optimal object illumination and can be dimmed.

1. Connect the digital microscope to the PC.
2. The brightness of the lighting can be adjusted by pressing the dimmer button  several times.

---

## 14 Connecting to a PC

Use of this digital microscope requires a connection via USB to a PC running the Windows operating system (Windows 7 or above) or macOS (10.13 or above). Follow the steps below for proper connection:

1. Download and install the HiView\* software on the PC.
2. Mount the digital camera as described.
3. Insert the USB-A plug of the USB connection cable into a free USB port of the PC. If the connection is successful, the LED lights up. The digital microscope is ready for use.
4. Via the HiView\* software, the images from the digital microscope can be viewed and saved. Furthermore, the images can be labeled and measured.
5. In the software, click on "Device" and then on "Open". Connect to the digital microscope found there (click on "Open"). If you have started the software first and then connect the digital microscope, you must click on "Refresh" so that the digital microscope can be found.

*\*Info about downloading the software can be found on page 2 in this manual.*

## 15 Adjust stand

Depending on the size of the observation object or when changing the magnification setting, the position of the microscope head must be adjusted in height. To do this, proceed as follows.

1. Carefully loosen the locking device on the side.
2. Set the desired new alignment and lock the locking screw again.

---

## 16 Adjusting the zoom

The microscope optics have a zoom range from 50x to 1000x.

In the software the zoom range can be changed by setting the display ratio from 100% to another value, e.g. 50% or 200%.

**NOTICE! When changing the zoom range, the focus may also need to be adjusted. (see also "Focusing the object")**

## 17 Focusing the object

For optimal image sharpness, the observation object must be brought into the focus range of the digital camera. To do this, proceed as follows:

1. Set the required position in the holder for the digital microscope and correct it slightly afterwards if necessary.
2. Set the desired magnification setting in the software (e.g. Display Ratio 100%).
3. Now turn the zoom wheel until the image appears sharp on the monitor.

## 18 Recording

Name: <input type="text" value="HD camera"/>	Here you select your digital microscope and establish a connection.
 Photo	Take photo
 Video	Record video
Resolution: <input type="text" value="640x480"/>	Set resolution (e.g. to 1920x1080 for the best picture quality)
 Files	Opens the save folder specified under settings.

Choose Ruler/Line/Text appearance. Choose save folder path and photo format. Change Language.

The following tools are available for labeling and measuring the exposures:

## 19 Technical data

Zoom range:	50x to 1000x (Display Ratio 50% - 200%)
Working distance:	15mm (50x) to 40mm (1000x)
Illumination:	8 LEDs
Digital camera:	fixed (CMOS sensor)
Image format:	BMP, JPG
Image Resolution:	640x480, 1280x720, 1920x1080** pixels
Video format:	AVI (30fps)
PC connection:	USB-A*
Power supply:	5V via USB connection*
*Data transfer and power supply are carried out via the USB port on the PC at the same time	
**Interpolated	

## 20 CE declaration of conformity

**CE** A "Declaration of conformity" in accordance with the applicable directives and corresponding standards has been prepared by Bresser GmbH. The full text of the CE declaration of conformity is available at the following Internet address: [www.bresser.de/download/9652200/CE/9652200\\_CE.pdf](http://www.bresser.de/download/9652200/CE/9652200_CE.pdf)

## 21 UKCA declaration of conformity

**UK  
CA** A "Declaration of Conformity" in accordance with the applicable directives and relevant standards has been issued by Bresser GmbH. The full text of the UKCA Declaration of Con-

---

formity is available at the following web address:  
[www.bresser.de/download/9652200/  
UKCA/9652200\\_UKCA.pdf](http://www.bresser.de/download/9652200/UKCA/9652200_UKCA.pdf)

**Bresser UK Ltd.** • Suite 3G, Eden House, Enterprise Way,  
Edenbridge, Kent TN8 6HF, Great Britain

## 22 Disposal

Dispose of the packaging materials according to its type. Information on proper disposal can be obtained from the municipal waste disposal service provider or environmental agency.

Do not dispose of electronic devices in the household garbage!

According to the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment and its transposition into national law, used electrical equipment must be collected separately and recycled in an environmentally sound manner.

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**CE** „Vyhlásenie o zhode“ v súlade s platnými smernicami a príslušnými normami vyhotovila spoločnosť Bresser GmbH: