Mikroskop · Microscope ·

Duolux / BioDiscover / Biolux / Erudit /
Biorit

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EN Instruction manual

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Contents

1 Imprint (German)	12
2 Validity note	12
3 About this Instruction Manual	12
4 Intended Use	12
5 General safety instructions	13
6 Parts overview	15
7 Choice of location	16
8 Setting up power supply and turning on the device	16
9 Inserting slides	16
10 Observation	16
11 Cleaning and maintenance	17
12 Disposal	17
13 Warranty	18
14 EC Declaration of Conformity	18

1 Imprint (German)

Bresser GmbH

Gutenbergstr. 2
46414 Rhede
Germany

<http://www.bresser.de>

For any warranty claims or service enquiries, please refer to the information on "Warranty" and "Service" in this documentation. We apologize for any inconvenience caused by the fact that we cannot process enquiries or submissions sent directly to the manufacturer's address.

Errors and technical changes excepted.

© 2020 Bresser GmbH

All rights reserved.

The reproduction of this documentation - even in extracts - in any form (e.g. photocopy, print, etc.) as well as the use and distribution by means of electronic systems (e.g. image file, website, etc.) without the prior written permission of the manufacturer is prohibited.

The designations and brand names of the respective companies used in this documentation are generally protected by trade, trademark and/or patent law in Germany, the European Union and/or other countries.

2 Validity note

This documentation is valid for the products with the following article numbers:

5116200 8855000 8855001 8855002 8855003 8855004 8855100 9039000 9039001 5012000
5013000 5110000 5101000 micro-bio-beginner

Manual version: v012021a

Manual designation:

Manual_Biological-Microscopes-Beginners_de-en_BRESSER_v012021a

Always provide information when requesting service.

3 About this Instruction Manual

NOTICE

These operating instructions are to be considered a component of the device.

Read the safety instructions and the operating manual carefully before using this device.

Keep this instruction manual in a safe place for future reference. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

4 Intended Use

- This device is not intended for use by persons (including children) with impaired physical, sensory or mental abilities, or for lack of experience and/or knowledge, unless they are supervised by or have received instructions from a person responsible for their safety on how to use this device.
- This device is for personal use only.
- It was developed for the magnified display of things in nature.
- The device is for indoor use only.

5 General safety instructions

DANGER

Risk of an electric shock!

This device contains electronic parts that are powered by a power source (AC adapter and/or batteries). Improper use of this product may result in electric shock. Electric shock can cause serious or fatal injuries. It is therefore imperative that you observe the following safety information.

- Disconnect the power supply by pulling the mains plug when the appliance is not in use, in case of a longer interruption of operation and before any maintenance and cleaning work.
- Place your device so that it can be disconnected from the power supply at any time. The power outlet should always be near your appliance and should be easily accessible, as the plug of the power cord serves as a disconnect device from the mains supply.
- To disconnect the unit from the mains, always pull the mains plug and never pull the cable!
- Check this device, cables and connections for damage before use.
- Never attempt to operate a damaged device, or a device with damaged electrical parts! Damaged parts must be replaced immediately by an authorized service agent.
- Operate the device only in a completely dry environment and do not touch the device with wet or damp body parts.

DANGER

Danger of suffocation!

Improper use of this product may result in suffocation, especially for children. It is therefore imperative that you observe the following safety information.

- Keep packaging materials (plastic bags, rubber bands, etc.) away from children! There is a DANGER OF CHOKING [These pose a CHOKING HAZZARD]
- This product contains small parts that can be swallowed by children! There is a DANGER OF CHOKING [These pose a CHOKING HAZZARD]
- If small parts has been swallowed, consult a doctor immediately!

DANGER

Explosion hazard!

Improper use of this product may result in fire. It is essential that you observe the following safety information in order to avoid fires.

- Do not expose the device to high temperatures. Use only the power supply provided. Never short-circuit the device or throw it into a fire. Excessive heat and improper handling of this product can result in short-circuits, fires or even explosions.

CAUTION

Danger of injury!

This device contains components and/or accessories that can cause minor to severe injuries in case of any improper use. Therefore please read the safety instructions below to avoid any bodily injury.

- Tools with sharp edges and points are often used when working with this device. Because there is a risk of injury from such tools, store this device and all tools and accessories in a location that is out of the reach of children.

-
- Children must not have access to the included chemicals and liquids. Do not drink the chemicals. Wash hands thoroughly with running water after using the chemicals. In the event that the chemicals come into contact with your eyes or mouth, rinse thoroughly with water. If you are in pain after exposure, contact a doctor immediately and take the substances with which you came into contact with you.

**CAUTION****Fire hazard!**

Improper use of this product may result in fire. It is essential that you observe the following safety information in order to avoid fires.

-
- Never cover the ventilation slots or cooling fins of the device while using it or as long as it has not sufficiently cooled down!

**NOTICE****Danger of material damage!**

Improper handling may result in damage to the unit and/or accessories. Therefore, use the device only in accordance with the following safety information.

-
- Do not disassemble the device! In the event of a defect, please contact your dealer. They will contact the Service Center and can arrange the return of this device for repair if necessary.
 - Do not expose the device to high temperatures and protect it from water and high humidity.
 - Do not subject the device to excessive vibrations.
 - Only use accessories and spare parts for this device that comply with the technical specifications.

6 Parts overview

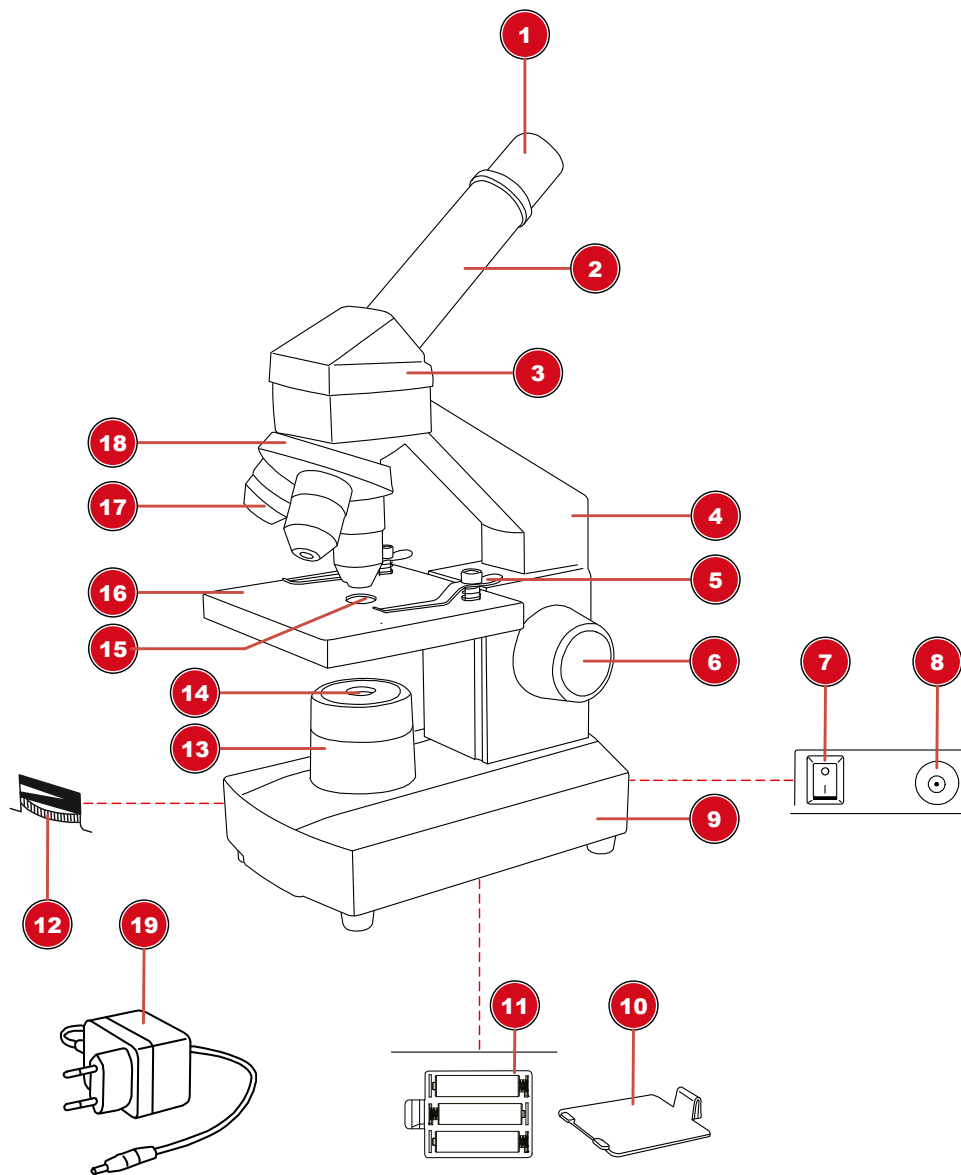
Illustration 2: Parts overview for biological microscopes

1 Eyepiece (can also be equipped with a pair of eyepieces) (optional equipment, depending on model)	2 Eyepiece holder
3 Microscope head	4 Microscope arm
5 Object fixing clamps	6 Focussing wheel
7 On/Off switch	8 DC socket for coaxial/barrel connector (can also be equipped with a C14 mains plug alternatively) (optional equipment, depending on model)
9 Microscope base	10 Battery compartment cover (optional equipment, depending on model)
11 Battery compartment (optional equipment, depending on model)	12 Rotary dimmer (optional equipment, depending on model)
13 Lighting unit	14 Lens
15 Light opening	16 Object stage
17 Objective	18 Objective nosepiece

19 Power adapter with Euro-plug, power cable and coaxial/barrel connector (can also be equipped with a Euro-plug and C13 mains socket) (optional equipment, depending on model)

7 Choice of location

Select a suitable location before installation and commissioning.

Place the unit on a stable, level and vibration-free surface.

8 Setting up power supply and turning on the device

Depending on the model, the power supply can be obtained by battery or mains power (power adapter/cable).

Battery operation (depending on model)

1. Remove the battery compartment cover.
2. Insert the batteries into the battery compartment.
3. Re-attach the battery compartment cover.
4. Move ON/OFF switch to position 'I' to turn on the device.
5. Turn the rotary dimmer to set the illumination brightness. (optional function, depending on model)
6. Move On/Off switch to 'O' position to turn off the device.

Mains operation (depending on model)

7. Put the DC connector of the power cable into the DC connection socket of the device.
8. Put the Euro-plug of the power cable into the wall socket (230V maximum voltage).
9. Move ON/OFF switch to position 'I' to turn on the device.
10. Turn the rotary dimmer to set the illumination brightness. (optional function, depending on model)
11. Move On/Off switch to 'O' position to turn off the device.

9 Inserting slides

1. Move the object plate downwards completely by turning the focus wheel.
2. Push down both object clamps at the backside that they lift up at the frontside.
3. Slide the microscope slide from the frontside below the raised object clamps.
4. Let the object clamps fall back carefully until the slide is fixed under them.

10 Observation

NOTICE

Using eyepieces and objectives

When using a specific eyepiece or objective, you will get the individual magnification of an object being observed. The magnification is imprinted on the eyepiece and objective with a large number (e.g. 10x for 10-fold magnification, 4 for 4-fold magnification).

1. Move the object plate downwards completely by turning the focus wheel.

2. Insert a slide and place it in the middle under the objective.
3. Set an objective with lowest magnification below the object plate.
4. Insert an eyepiece with lowest magnification in the eyepiece holder.
5. Turn on the device and the illumination.
6. Look through the eyepiece and move the object plate upwards by turning the focus wheel until a sharp image is visible.
7. If necessary, adjust the image sharpness by turning the fine focus adjustment.

NOTICE

Avoid damages to the specimen!

Make sure you turn the focus wheel very slowly, to avoid the objective touch the slide or specimen, and thus to avoid damages to either of them, or to both. Only turn the focus wheel until it stops and do not apply excessive pressure.

11 Cleaning and maintenance

- Protect the device from dust and moisture!
- Avoid fingerprints and similar contamination of the optical surfaces.
- The residual moisture must be completely degraded before use.
- Disconnect the power supply by pulling the mains plug when the appliance is not in use, in case of a longer interruption of operation and before any maintenance and cleaning work.
- Only clean the device externally using a dry cloth. Do not use cleaning solution to prevent damage to the electronic parts.
- Do not use organic solvents (e.g. alcohol, ether, acetone, xylol or other dilutions) to clean lacquered parts or plastic parts!
- Disassemble optical elements (e.g. objective, eyepiece, etc) before cleaning.
- Blow away loose dust from the lens surfaces first.
- Use high quality lens cleaning towels or soft cloth and moisten it with a little bit of pure alcohol (available in a drug store). Wipe the lens surface with it.
- Pull a dust protection cover* over the microscope and store it in a closed container at a dry and mould-free place. Store the microscope and the accessories in the relevant containers when they are not used for a longer time. It is recommended to store also objectives and eyepieces in closed containers with desiccant.

(*depending on the model a dust protection cover and box can be part of the content)

NOTICE

Remember:

A well maintained microscope will keep its optical quality for years and thus maintain its value.

12 Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

According to the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment and its transposition into national law, used electrical equipment must be collected separately and recycled in an environmentally sound manner.

Batteries and rechargeable batteries must not be disposed of with household waste. You are legally obliged to return used batteries and accumulators and can return the batteries after use either in our sales outlet or in the immediate vicinity (e.g. in the trade or in municipal collection points) free of charge.

Batteries and accumulators are marked with a crossed-out dustbin and the chemical symbol of the pollutant, "Cd" stands for cadmium, "Hg" stands for mercury and "Pb" stands for lead.

Cd¹

Hg²

Pb³

13 Warranty

The regular guarantee period is 5 years and begins on the day of purchase. You can consult the full guarantee terms and details of our services at www.bresser.de/warranty_terms.

14 EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. This can be viewed any time upon request. <http://www.bresser.de> • info@bresser.de

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

e-mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd.
Suite G3, Eden House
Enterprise Way
Edenbridge, Kent TN8 6HF
United Kingdom

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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Folux B.V.